



			SacI				XhoI				HindIII	EcoRI			Sall	
1315	<u>End mCherry</u>	<u>BspEI</u>	<u>CTC</u>	<u>BglII</u>	<u>AGA TCT</u>	<u>CGA GCT</u>	<u>CAA GCT</u>	<u>TCG AAT</u>	<u>TCT GCA</u>	<u>GTC</u>						
		<u>TAC AAG</u>	<u>TCC GGA</u>													
	<u>Sall</u>	<u>KpnI</u>	<u>SacII</u>	<u>XmaI</u>												
1363	<u>AccI</u>	<u>Asp718</u>	<u>GCG GGC</u>	<u>CGG GGA</u>	<u>TCC ACC</u>	<u>GGA TCC</u>	<u>ACC GGA</u>	<u>TCT AGA</u>	<u>TAA CTG</u>	<u>ATC ATA</u>						
	<u>GAC</u>	<u>GGT ACC</u>		<u>Apal</u>	<u>BamHI</u>			<u>XbaI*</u>		<u>BclI</u>						

\* The XbaI site is methylated in the DNA provided by Clontech. If you wish to digest the vector with XbaI enzyme, you will need to transform the vector into a dam- host and make fresh DNA.

#### pmCherry-C1 Restriction Map and Multiple Cloning Site (MCS).

#### Description

pmCherry-C1 is a mammalian expression vector designed to express a protein of interest fused to the C-terminus of mCherry, a mutant fluorescent protein derived from the tetrameric *Discosoma sp.* red fluorescent protein, DsRed (1). The excitation and emission maxima of the native mCherry protein are 587 nm and 610 nm, respectively. Expression of fusion proteins that retain the fluorescent properties of the unmodified mCherry protein can be monitored by flow cytometry and their localization *in vivo* can be determined by fluorescence microscopy.

The multiple cloning site (MCS) in pmCherry-C1 is positioned downstream of the mCherry coding sequence. A Kozak consensus sequence is located immediately upstream of the mCherry gene to enhance translational efficiency in eukaryotic systems (2). SV40 polyadenylation signals downstream of the mCherry gene and the MCS direct proper processing of the 3' end of the mCherry (or fusion gene) mRNA.

The vector backbone contains an SV40 origin for replication in mammalian cells expressing the SV40 large T antigen, a pUC origin of replication for propagation in *E. coli*, and an f1 origin for single-stranded DNA production. A neomycin-resistance cassette (Neo<sup>r</sup>) allows stably transfected eukaryotic cells to be selected using G418. This cassette consists of the SV40 early promoter ( $P_{SV40\ e}$ ), the Tn5 neomycin/kanamycin resistance gene, and polyadenylation signals from the herpes simplex virus thymidine kinase (HSV TK) gene. A bacterial promoter ( $P_{Kan^r}$ ) upstream of the cassette confers kanamycin resistance in *E.*

(PR8X2616; published 24 October 2008)



**Clontech**

United States/Canada  
800.662.2566

Asia Pacific  
+1.650.919.7300

Europe  
+33.(0)1.3904.6880

Japan  
+81.(0)77.543.6116

Clontech Laboratories, Inc.  
A Takara Bio Company  
1290 Terra Bella Ave.  
Mountain View, CA 94043  
Technical Support (US)  
E-mail: tech@clontech.com  
www.clontech.com

*coli*.

### Use

The gene of interest must be cloned into pmCherry-C1 so that it is in-frame with the mCherry coding sequence. The gene must include an initiation codon (ATG), and lack in-frame stop codons.

The pmCherry-C1 vector can be transfected into mammalian cells using any standard transfection method. If required, stable transfectants can be selected using G418 (3). pmCherry-C1 can also be used as a cotransfection marker, as the unmodified vector will express mCherry in mammalian cells.

For Western analysis, either the Living Colors® DsRed Polyclonal Antibody (Cat. No. 632496) or the DsRed Monoclonal Antibody (Cat. Nos. 632392 and 632393) can be used to detect the mCherry protein.

### Location of features

- $P_{CMVIE}$  (human cytomegalovirus immediate early promoter): 1–589  
Enhancer region: 59–465; TATA box: 554–560  
Transcription start point: 583  
C→G mutation to remove the SacI site: 569
- mCherry (human codon-optimized)  
Kozak consensus translation initiation site: 606–616  
Start codon (ATG): 613–615; Stop codon: 1395–1397, 1399–1401 & 1403–1405  
Last amino acid: 1318–1320
- MCS (multiple cloning site): 1321–1408
- SV40 early polyA<sup>+</sup> signals  
Polyadenylation signals: 1541–1546 & 1570–1575; mRNA 3' ends: 1579 & 1591
- f1 origin of replication: 1638–2093 (complementary)
- $P_{Kan^r}$  (bacterial promoter for Kan<sup>r</sup> gene expression)  
–35 region: 2155–2164; –10 region: 2178–2183  
Transcription start point: 2190
- $P_{SV40e}$  (SV40 early promoter and enhancer sequences)  
Enhancer (72-bp tandem repeats): 2267–2338 & 2339–2410  
21-bp repeats: 2414–2434, 2435–2455 & 2457–2477  
Early promoter element: 2490–2496  
Major transcription start points: 2486, 2524, 2530 & 2535
- SV40 origin of replication: 2434–2569
- Kan<sup>r</sup>/Neo<sup>r</sup> (kanamycin/neomycin resistance gene)  
Neomycin phosphotransferase coding sequences:  
Start codon (ATG): 2618–2620; stop codon: 3410–3412  
G→A mutation to remove the PstI site: 2800  
C→A (Arg to Ser) mutation to remove BssHIII site: 3146
- HSVTK polyA<sup>+</sup> (herpes simplex virus thymidine kinase polyadenylation signals): 3648–3653 & 3661–3666
- pUC origin of replication: 3997–4640

### Propagation in *E. coli*

- Suitable host strains: DH5 $\alpha$ , HB101, and other general purpose strains. Single-stranded DNA production requires a host containing an F plasmid such as JM109 or XL1-Blue.
- Selectable marker: plasmid confers resistance to kanamycin (50  $\mu$ g/ml) in *E. coli* hosts.
- *E. coli* replication origin: pUC
- Copy number: high
- Plasmid incompatibility group: pMB1/Col E1

### Excitation and emission maxima of mCherry

- Excitation maximum = 587 nm
- Emission maximum = 610 nm

### References

1. Shaner, N. C., *et al.* (2004) *Nature Biotech.* **22**(12):1567-72.
2. Kozak, M. (1987) *Nucleic Acids Res.* **15**(20):8125–8148.
3. Gorman, C. (1985) In *DNA Cloning: A Practical Approach, Vol. II*. Ed. D. M. Glover (IRL Press, Oxford, U.K.) pp. 143–190.

**Note:** The attached sequence file has been compiled from information in the sequence databases, published literature, and other sources, together with partial sequences obtained by Clontech Laboratories, Inc. This vector has not been completely sequenced.

**Notice to Purchaser**

Clontech products are to be used for research purposes only. They may not be used for any other purpose, including, but not limited to, use in drugs, *in vitro* diagnostic purposes, therapeutics, or in humans. Clontech products may not be transferred to third parties, resold, modified for resale, or used to manufacture commercial products or to provide a service to third parties without written approval of Clontech Laboratories, Inc.

The CMV promoter is covered under U.S. Patent Nos. 5,168,062, and 5,385,839 assigned to the University of Iowa Research Foundation.

DsRed-Express: Patent Pending.

Fruit Fluorescent Protein Products pmCherry, pmRaspberry, pmPlum, pmBanana, pmOrange, pmStrawberry and their variants:

Not-For-Profit Entities: Orders may be placed in the normal manner by contacting your local representative or Clontech Customer Service at 650.919.7300. At its discretion, Clontech grants Not-For-Profit Entities a non-exclusive, personal, limited license to use this product for non-commercial life science research use only. Such license specifically excludes the right to sell or otherwise transfer this product, its components or derivatives thereof to third parties. No modifications to the protein coding sequence may be made without express written permission from Clontech. Any other use of this product requires a license from Clontech. For license information, please contact a licensing representative by phone at 650.919.7320 or by e-mail at [licensing@clontech.com](mailto:licensing@clontech.com).

For-Profit Entities wishing to use this product are required to obtain a license from Clontech. For license information, please contact a licensing representative by phone at 650.919.7320 or by e-mail at [licensing@clontech.com](mailto:licensing@clontech.com).

This product is the subject of U.S. patents.

Clontech, the Clontech logo and all other trademarks are the property of Clontech Laboratories, Inc., unless noted otherwise. Clontech is a Takara Bio Company. ©2008 Clontech Laboratories, Inc.